

DASYLab and instruNet devices - more than 8 instruNet channels on worksheet will not save correctly

If you are working with DASYLab V13 or earlier and have more than 8 instruNet channels, the channel settings might not save properly in the DASYLab .dsb worksheet file.

To fix this, you will need to download and copy a new "Inet.dll" # v13.0.1.183 to the installed DASYLab file location. On Windows 64-bit systems it is ..\Program Files (x86)\DasyLab 13.0\. On 32-bit Windows systems it is ..\Program Files\DASYLab 13.0\ . Note that earlier DASYLab versions will have the version number in the folder name.

You will need Administrator privileges on Windows 7 or later to replace the file.

Note: This will replace the the DASYLab driver DLL, not the instruNet system drivers (iNet32.dll).

To download this file, visit:

http://gwinst.com.previewc28.carrierzone.com/ftp/DASYLab/inet_dll/v13-0-1-183/

It is documented at:

<http://www.gwinst.com/software/dasylab/installation.html#newdll>

Measurement Computing Data Acquisition Knowledgebase
<https://kb.mccdaq.com/KnowledgebaseArticle50598.aspx>